

Zone samplers | for all kinds of bulk goods



1. Multi sampler (closed chambers);
2. All layer sampler (open inner tube);
3. Uno sampler (one closed chamber).

Bürklee zone samplers are used for taking samples of bulk goods of any kind.

The big advantage of zone samplers is that all-layer samples can be collected from all layers of the container.

Various bulk goods can be collected with the lances, from the finest powder to large grains such as corn or nuts.

The zone samplers are available in four lengths. 55 cm (especially suitable for sacks), 85 cm long (especially for barrels, tubs, mixed containers) and 150 cm as well as 250 cm for big bags, silos, tanks.

Safe and reliable sampling is made even easier with useful special accessories such as sealing sleeves, Quali-Paper, cleaning brushes, emptying devices and close-it control seals.

The zone samplers are available in different models for different sampling tasks.

- ▶ Smallest possible gap between the inner and the outer tube
- ▶ Easy cleaning, removable tip
- ▶ Colour coding (Multi and Uno sampler)
- ▶ Ultra-pure materials
- ▶ Excellent processing, robust model



Colour coding: Identification to prevent mixing up multi samplers and single samplers.



contact@multilab.ro service@multilab.ro info@multilab.ro

Hot-Line 0724044800

Rel.Client 0730 711101

Desfacere 072 8930531 / 073 5876690 / 0735876689 / 0728930530

Service 0728930535 / 0728930536 / 0731309088

Secretariat 021 3127144 / 45 Fax. 021 3127141

Zone samplers | for all kinds of bulk goods



All-layer sampler, with open inner tube, no separate chambers, and large sampling volume. For all-layer samples through all the layers of the sample.

- ▶ Aluminium or V4A (1.4404)
- ▶ 25 mm Ø

Corresponds to ISTA standard for seed lengths of 10 mm (wheat, oats) and 5 mm (flax, vetches).



Multi sampler with several closed chambers, for multi point sampling from several predetermined sampling depths.

- ▶ Aluminium anodised, V4A (1.4404) or V4A/PTFE
- ▶ Version made of V4A/PTFE with low weight due to PTFE sample chambers.
- ▶ With colour coding system
- ▶ 25 mm Ø



Uno sampler made of stainless steel V4A/PTFE with a closed sample chamber for point sampling from a predetermined sampling depth.

- ▶ Outer tube stainless steel V4A (1.4404), sample chambers made of PTFE
- ▶ With colour coding system
- ▶ 25 mm Ø



close-it seals openings. See page 88.

All-layer-sampler

Material	Total length cm	Insertion depth mm	No. of openings	Chamber volume ml	Price €	Item no.
aluminium	55	430	3	70		5317-2055 N4
aluminium	85	710	3	130		5317-2085 N4
aluminium	150	1355	5	260		5317-2150 N4
V4A	55	430	3	70		5316-2055 N4
V4A	85	710	3	130		5316-2085 N4
V4A	150	1355	5	260		5316-2150 N4

Multi sampler

Material	Total length cm	Insertion depth mm	Number of collection chambers	Chamber volume ml	Price €	Item no.
aluminium	55	430	3	14		5317-1058 N4
aluminium	85	710	3	17		5317-1085 N4
aluminium	150	1355	5	17		5317-1150 N4
V4A	55	430	3	14		5316-3055 N4
V4A	85	710	3	17		5316-3085 N4
V4A	150	1355	5	17		5316-3150 N4
V4A/PTFE	55	430	3	14		5316-0055 N4
V4A/PTFE	85	710	3	17		5316-0085 N4
V4A/PTFE	150	1355	5	17		5316-0150 N4

Uno sampler

Material	Total length cm	Insertion depth mm	Number of collection chambers	Chamber volume ml	Price €	Item no.
V4A/PTFE	55	430	1	17		5316-1055 N4
V4A/PTFE	85	710	1	17		5316-1085 N4
V4A/PTFE	150	1355	1	17		5316-1150 N4
V4A	150	1355	1	17		5316-4150 N4

Use

1. Turn the handle, close the chambers.
2. Insert the sampler to the required depth.
3. Open the chambers, sample flows into the chamber.
4. Close the chambers.
5. Pull out the sampler.
6. Close the insertion opening with the close-it control seal.
7. Empty the chambers onto Quali-Paper or with the emptying device.

Zone samplers | for all kinds of bulk goods



Jumbo sampler with a tube diameter of 50 mm, ideal for coarse grains and larger quantities.

With an open inner tube and large sampling volume, for all-layer samples through all layers of the sample or with sample chambers separated by PTFE separators for multi point sampling from several pre-determined sampling depths.

- ▶ Aluminium or aluminium/PTFE
- ▶ 50 mm Ø

Jumbo sampler

Model	Total length cm	Material	Insertion depth mm	No. of openings	Chamber volume ml	Price €	Item no.
All layer - open inner tube	85	aluminium	710	3	880		5319-1085 N4
All layer - open inner tube	150	aluminium	1355	5	1700		5319-1150 N4
All layer - open inner tube	250	aluminium	2355	7	2900		5319-1250 N4
Multi - closed chambers	85	aluminium/PTFE	710	3	390		5319-2085 N4
Multi - closed chambers	150	aluminium/PTFE	1355	5	650		5319-2150 N4
Multi - closed chambers	250	aluminium/PTFE	2355	7	910		5319-2250 N4



Accessories for zone samplers



Emptying device



Zone sampler sealing sleeves



Quali-Paper for zone sampler



Cleaning brushes

Emptying device

Model	Material	Length cm	Price €	Quantity scale from	Scale price €	Item no.
device with 3 funnels	V2A	100		5		5317-0030 N4
emptying funnel separate	V2A			3		5317-0033 N4

Sealing sleeves

Dia. mm	Pieces per pack	Price €	Quantity scale from	Scale price €	Item no.
25	3		10		5317-0011 N4
50	3		10		5317-0010 N4

Quali-Paper

L x W cm	Pcs. per pack	Price €	Quantity scale from	Scale price €	Item no.
100 x 23	50		10		5317-0001 N4

Cleaning brushes

Dia. mm	Length cm	Material	Price €	Quantity scale from	Scale price €	Item no.
25	40	PVC/stainless steel		10		5304-0040 N4
25	60	PVC/stainless steel		10		5304-0060 N4
25	80	PVC/stainless steel		10		5304-0080 N4
25	100	PVC/stainless steel		10		5304-0100 N4
50	100	PVC/stainless steel		5		5304-5100 N4
50	300	PVC/stainless steel		5		5304-5300 N4

Zone sampler Spiralus



- ▶ Representative sample of the container layers.
- ▶ Aluminium, anodised
- ▶ 38 mm Ø
- ▶ Low weight

Guarantees precise zone sampling through all layers of the sample. The spiral arrangement of the sampling slits means that these are opened one after the other from bottom to top. In this way the lowest opening is filled first and then those above it in sequence until the complete Spiralus has been filled.

The shape of the sampling slits prevents seeds from being cut up. The Spiralus provides an excellent all-layer sample.

Applications


Spiralus is particularly suitable for sampling coarse-grained bulk goods, seeds and pellets at locations such as soil heaps, warehouses, trucks, railway goods wagons, etc.

Zone sampler Spiralus					
Total length cm	No. of openings	Weight g	Chamber volume ml	Price €	Item no.
145	8	1400	770		5318-0145 N4
200	8	2000	1100		5318-0200 N4
Accessories					
Description	Price €	Item no.			
Cleaning brushes 50 mm dia., length 1000 mm		5304-5100 N4			

EasySampler | easy to empty



- ▶ Stainless steel V4A (1.4404)
- ▶ For sample opening from 52 mm Ø
- ▶ Length of the sample chamber 500 mm

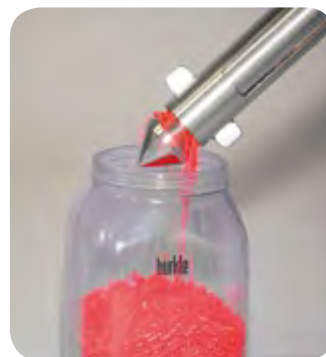

It is possible to take multiple samples from bulk goods with this sampler quickly and easily.

Emptying has been greatly simplified and in this model takes place via the tip. The advantage: The sample can be emptied directly into the sample container without having to disassemble the EasySampler, in order to obtain a mixed sample for instance.

The bulk material must be free-flowing or self-flowing.

EasySampler has been developed in cooperation with an international chemical concern based in Ludwigshafen, Germany.

EasySampler				
Tube Ø mm	Total length cm	Volume ml	Price €	Item no.
32	104,5	200		5360-1000 N2

All-layer-scratcher



The scratcher is a special sampler for powders and granules that do not flow freely. The sample is transferred to the closable inner tube using a scraper. Sliding PTFE separators are used to divide the inner tube into different zones. The separators also prevent the sample from being mixed.

- For caked bulked goods
- Zones can be varied due to moveable PTFE separators.
- Made of stainless steel V4A (1.4404) and PTFE

All-layer-scratcher			
			  
Dia. mm	Length mm	Price €	Item no.
25	1000		5351-1000 N2
 Accessories			
Description	Price €	Item no.	
Cleaning brushes 25 mm dia., length 100 cm		5304-0100 N4	



contact@multilab.ro service@multilab.ro info@multilab.ro

Hot-Line 0724044800

Rel.Client 0730 711101

Desfacere 072 8930531 / 073 5876690 / 0735876689 / 0728930530

Service 0728930535 / 0728930536 / 0731309088

Secretariat 021 3127144 / 45 Fax: 021 3127141

Scratcher | for non-pourable substances






The scratcher has particularly been designed for sampling cohesive, caked and not free-flowing substances.

- Stainless steel V4A (1.4404)
- Electro-polished
- Total length 1000 mm

Use

1. Insert the closed lance into the sample.
2. Open the inner tube by turning the sample chamber when at the required depth.
3. Turn the scratcher by 360° in the sample, the scraper directs the caked material into the sample chamber.
4. Close the sample chamber.

Scratcher			
			  
Chamber volume ml	Flute length mm	Outer dia. with scraper mm	Price €
5	22	25	
10	45	25	
25	110	25	
			Item no.
			5347-0005 N2
			5347-0010 N2
			5347-0025 N2

Quality assurance starts with the inspection of incoming goods! An all-layer sample of bulk goods is no longer a problem with PowderProof®, the sampler for powders and granules.


PowderProof® is the practically proven professional solution for mobile sampling of powders and granules. Different lances and augers can be attached to the robust drive unit and swapped without any problems. This modular design makes PowderProof® a versatile instrument for sampling powders and granules from drums, barrels, bags or sacks.

Safe and reliable sampling is made even easier with useful accessories such as basket-enclosed bottles, sample bags and close-it control seals.

PowderProof® with lance for horizontal use, such as sampling from sacks, with lance for vertical use, for sampling from open containers such as barrels, drums or bags.




PowderProof® | mobile powder sampling

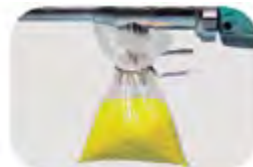

- Many uses
- Stainless steel V4A (1.4404)
- Incl. motor 12 V, two re-chargeable batteries, fast charger, clip for holding the bag and cleaning brush
- Additional augers available as accessory

Use

1. The lance is inserted into the sack/container with the spiral screw rotating.
2. The powder or granules are immediately transported through the lance and filled directly into a sample bag or bottle at the lance outlet.
3. Any excess sample remaining in the lance is returned to the sack by reversing the spiral screw direction

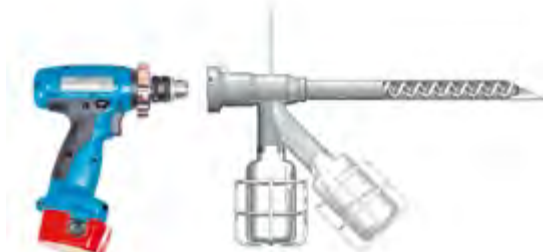

PowderProof for horizontal and vertical use

PowderProof®

Model	Lance	Length cm	Price €	Item no.
for powders up to 2mm dia.	horizontal	30		5302-0302 N2
for powders up to 2mm dia.	vertical	60		5302-0602 N2
for powders up to 2mm dia.	vertical	90		5302-0902 N2
for powders up to 2mm dia.	vertical	120		5302-1202 N2
for granules from 2 up to 4mm dia.	horizontal	30		5302-0301 N2
for granules from 2 up to 4mm dia.	vertical	60		5302-0601 N2
for granules from 2 up to 4mm dia.	vertical	90		5302-0901 N2
for granules from 2 up to 4mm dia.	vertical	120		5302-1201 N2

Accessories

Description	Price €	Item no.
Spiral screw for PowderProof for powders up to 2mm dia., 30 cm		5302-0312 N2
Spiral screw for PowderProof for powders up to 2mm dia., 60 cm		5302-0612 N2
Spiral screw for PowderProof for powders up to 2mm dia., 90 cm		5302-0912 N2
Spiral screw for PowderProof for powders up to 2mm dia., 120 cm		5302-1212 N2
Spiral screw for PowderProof for granules from 2 up to 4mm dia., 30 cm		5302-0311 N2
Spiral screw for PowderProof for granules from 2 up to 4mm dia., 60 cm		5302-0611 N2
Spiral screw for PowderProof for granules from 2 up to 4mm dia., 90 cm		5302-0911 N2
Spiral screw for PowderProof for granules from 2 up to 4mm dia., 120 cm		5302-1211 N2
Bag clip for PowderProof		0006-6259 N4
Transport case L x W 121 x 22,5 cm		3600-2120 N4
Sampling bottle LDPE 500 ml		0318-0500 N3
Basket for sample bottles		5302-2050 N2
Packaging bags with closure, 170x120 mm, 400 ml, 100 pieces		2348-0005 N3

Novartos | pharma zone sampler according to FDA guidelines for unit-dose sampling




The Novartos Multi has 14 sample chambers for up to seven sample zones for an all-layer sample.

Novartos was developed in co-operation with a globally-operating pharmaceuticals and chemicals corporation located in Basel, Switzerland.

- ▶ The sample inserts fit precisely and can easily be replaced.
- ▶ Inserts are available in volumes from 0 to 4.0 ml
- ▶ The insertion depths can easily be read from the laser-engraved depth scale
- ▶ Stainless steel V4A (1.4404/1.4571)
- ▶ Groove-free and hand-polished
- ▶ 25 mm Ø



Novartos Uno sampler with two sample chambers, for one sample zone.

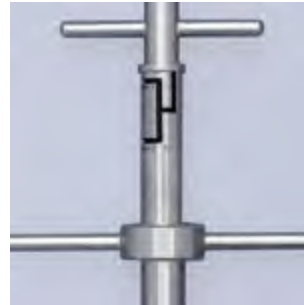

Novartos double-sample



Computer optimized inserts



Please order inserts separately!

Novartos the pharma zone sampler according to FDA Guidelines Unit-Dose-Sampling. It is now possible to withdraw a double-sample without inserting the lance repeatedly into the bulk material. Mixing is mainly prevented in this way.

One significant advantage of Novartos is the easy cleaning.

Complete filling and emptying of sample chambers has been improved substantially using computer-optimised geometry. There are no crevices or dead spaces where sampling residues can collect.

- Novartos Uno for targeted sampling
- Novartos Multi enables all-layer samples

Accessories

Description	Price €	Item no.
Inserts for Novartos Multi, capacity 0,0 ml		5358-0000 N4
Inserts for Novartos Multi, capacity 0,5 ml		5358-0005 N4
Inserts for Novartos Multi, capacity 1 ml		5358-0010 N4
Inserts for Novartos Multi, capacity 1,5 ml		5358-0015 N4
Inserts for Novartos Multi, capacity 2 ml		5358-0020 N4
Inserts for Novartos Uno, capacity 0 ml		5358-0100 N4
Inserts for Novartos Uno, capacity 0,5 ml		5358-0105 N4
Inserts for Novartos Uno, capacity 1 ml		5358-0110 N4
Inserts for Novartos Uno, capacity 2 ml		5358-0120 N4
Inserts for Novartos Uno, capacity 3 ml		5358-0130 N4
Inserts for Novartos Uno, capacity 4 ml		5358-0140 N4

Novartos Multi

Model	Length mm	Price €	Item no.
without inserts	1000		5358-1000 N4

Accessories

Description	Price €	Item no.
Transport case L x W 121 x 22,5 cm		3600-2120 N4

Novartos Uno

Model	Length mm	Price €	Item no.
without inserts	600		5358-0600 N4
without inserts	1000		5358-1001 N4
without inserts	1400		5358-1400 N4

Accessories

Description	Price €	Item no.
Transport case L x W 121 x 22,5 cm		3600-2120 N4

MicroSampler | useful for taking small samples, sterilisable



Use

1. Insert MicroSampler into powder.
2. At the predetermined point, open the volume tip by pressing the button, the sample flows in.
3. Close the volume tip, the sample is collected.
4. Pull back the lance, open the volume tip again, empty the sample.

The MicroSampler has especially been designed for the sampling of small quantities of powder.


Volumes of 0.2 ml, 0.5 ml, 1.0 ml, 2.0 ml, 5.0 ml and 10 ml can be sampled easily, quickly and cleanly from different sampling depths from sacks or containers.

The tips of the MicroSampler are interchangeable. This means that tips with different volumes can be used.

- Point sampling
- Various lengths and volumes are available
- Tip can be removed, easy cleaning
- Stainless steel V4A (1.4404)
- Sterilisable
- Additional tips are available in different sizes



Please order lance and head separately!

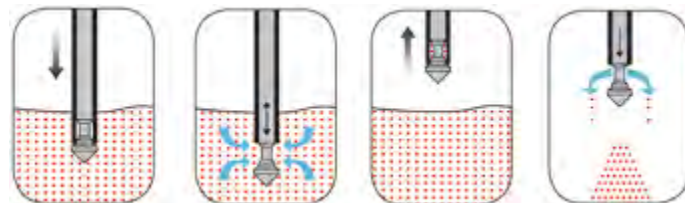

MicroSampler




Tube Ø mm	Length cm	Model	Price €	Item no.
12	55	without head		5307-1055 N4
12	85	without head		5307-1085 N4
12	120	without head		5307-1120 N4
25	55	without head		5307-2055 N4
25	85	without head		5307-2085 N4
25	120	without head		5307-2120 N4

Accessories

Description	Price €	Item no.
MicroSampler heads fits tube with 12 mm dia., capacity 0,2 ml		5307-0102 N4
MicroSampler heads fits tube with 12 mm dia., capacity 0,5 ml		5307-0105 N4
MicroSampler heads fits tube with 12 mm dia., capacity 1ml		5307-0110 N4
MicroSampler heads fits tube with 25 mm dia., capacity 2 ml		5307-0202 N4
MicroSampler heads fits tube with 25 mm dia., capacity 5 ml		5307-0205 N4
MicroSampler heads fits tube with 25 mm dia., capacity 10 ml		5307-0210 N4

PharmaPicker | for sampling very small quantities

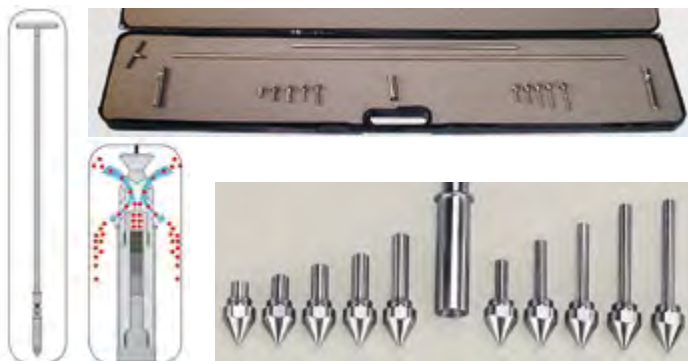

For sampling very small quantities of highly potent, expensive or active powders for analysis. The smallest sample quantities from 0.1 ml to 2.5 ml can be collected with the PharmaPicker. By simply exchanging the volume tip in the collection cylinder for 0.1 - 1.2 ml or 1.25 - 2.5 ml, the sampling volume is modified.

The PharmaPicker has a length of 60 cm and can be extended repeatedly by 50 cm or 100 cm, maximum length 3.50 m. The PharmaPicker is available separately or as PharmaPicker Case in a handy transport box.

PharmaPicker Case consisting of: Collection cylinders in both sizes, extension rod 50 cm and 100 cm, handle, volume tips 0.4 ml, 0.6 ml, 0.8 ml, 1.0 ml, 1.2 ml, 1.4 ml, 1.8 ml, 2.0 ml, 2.3 ml and 2.5 ml (alternative sample volumes on request), transport box.

- ▶ Stainless steel V4A
- ▶ electro-polished

Please order lance and head separately!

PharmaPicker						
Model	Price €	Item no.				
with collection cylinder for 0.10 - 1.20 ml, without head		5346-1000 N2				
with collection cylinder for 1.25 - 2.50 ml, without head		5346-2000 N2				
Case, completely		5346-0010 N2				

Accessories

Description	Price €	Item no.
Volume heads for PharmaPicker, capacity 0,10 ml		5346-1010 N2
Volume heads for PharmaPicker, capacity 0,20 ml		5346-1020 N2
Volume heads for PharmaPicker, capacity 0,60 ml		5346-1060 N2
Volume heads for PharmaPicker, capacity 1,00 ml		5346-1100 N2
Volume heads for PharmaPicker, capacity 2,00 ml		5346-2200 N2
Volume heads for PharmaPicker, capacity 2,50 ml		5346-2250 N2


Accessories

Description	Price €	Item no.
Extension rod PharmaPicker 500 mm		5380-0506 N4
Extension rod PharmaPicker 1000 mm		5380-1006 N4
Collection cylinder for 0.10 - 1.20 ml, without volume heads		5346-1001 N2
Collection cylinder for 1.25 - 2.50, without volume heads		5346-2001 N2
Handle PharmaPicker		0006-8113 N3

Use

1. Insert the required volume tip into the collection cylinder.
2. Insert the PharmaPicker, the disk sleeve closes the collection cylinder automatically.
3. The disk sleeve opens the collection cylinder at the required point, and the desired sample is collected.
4. Withdraw the PharmaPicker completely.
5. Empty the sample directly from the collection cylinder into the analysis container.

PharmaSpoon | for ultrapure sampling

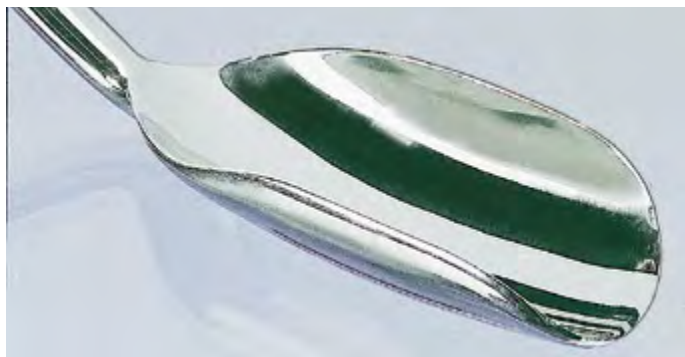

PharmaSpoon for taking samples of powdery substances, intermediate products and material in the pharmaceutical industry. Special materials with excellent surface qualities must be used for this to prevent cross-contamination. Our PharmaSpoon fulfils these requirements. Made of stainless steel V4A (1.4404), surface highly polished by hand (not electro-polished!). The excellent surface properties of the PharmaSpoon provide significant advantages.

- ▶ Possible residues are even visible in the trace sector (μg sector)
- ▶ The surface is easy to clean
- ▶ The PharmaSpoon is also suitable for sterile sampling



The PharmaSpoon was developed by an internationally known pharmaceutical company in Basel, Switzerland, and its usefulness has been confirmed in actual practice.

PharmaSpoon

Capacity ml	Total length mm	Price €	Quantity scale from	Scale price €	Item no.	
500	650		5		5324-0001 N2	

StickProof



Sampler for powder in the pharma area, made of stainless steel V4A, material 1.4404. The especially slender tip can be inserted easily and deeply in order to take the sample. The conical inlet of the tip allows the collection of either very small or larger quantities. StickProof is constructed in one piece without edges, grooves or crevices for perfect and easy cleaning. The sample can be filled directly into the sampling bag from the outlet end. Fasten the bag with a sturdy clamp.



Corresponds to ISTA standard for seed lengths of 10 mm (wheat, oats) and 5 mm (flax, vetches).

- ▶ Surface electro-polished
- ▶ Various sample quantities up to approx. 50 ml
- ▶ Incl. clamp and 100 PE sampling bags 120 x 170 mm

StickProof

Length mm	Dia. mm	Price €	Item no.	
410	25		5338-0410 N4	
1100	25		5338-1100 N4	

Accessories

Description	Price €	Item no.
StickProof clamp		5338-0001
Packaging bags with closure, 170x120 mm, 400 ml, 100 pieces		2348-0005 N3
close-it control seals see page 88		

Sampling lance Milky | acc. to EN ISO 707



Universal powder and granules sampler made of stainless steel V4A. Developed initially for milk powder samples, conforms to standards of DIN and IDF (International Dairy Federation). The collector is inserted into bags or casks and the sample is removed using a semi-circular profile lance.



- ▶ Stainless steel V4A (1.4404)
- ▶ Insertion depth 385 mm
- ▶ Total length 480 mm
- ▶ Outer diameter approx. 28 mm


close-it and close-it food

See page 88.

Sampling lance Milky

Volume ml	Price €	Quantity scale from	Scale price €	Item no.
150		5		5321-0300 N4

Sampler Tubus

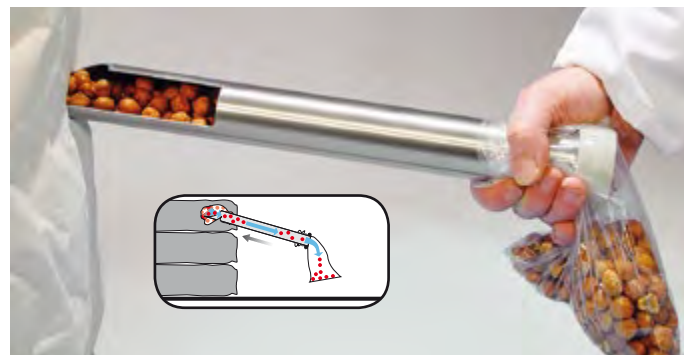

Sampler for pourable bulk goods up to approx. 1 cm in diameter. This makes it possible, for example, to remove hazel nuts, blends of tea, grain or similar items from large sacks.

- ▶ Also for coarse-grained bulk goods
- ▶ Made of stainless steel V4A (1.4404)
- ▶ Special geometry of the tip

Corresponds to ISTA standard for seed lengths of 10 mm (wheat, oats) and 5 mm (flax, vetches).

Use

1. Insert Tubus into a sack to the desired sampling depth, while keeping the collection tube closed using your hand.
2. When you remove your hand, the sample flows directly through the open tube into the container or sampling bag beneath it.


Sampler Tubus

Length cm	Dia. mm	Price €	Quantity scale from	Scale price €	Item no.
40	40		10		5382-0040 N4
80	40		10		5382-0080 N4

Accessories

Description	Price €	Item no.
close-it control seals see page 88		

QuickPicker | collects samples quickly and channels them directly into the sampling bottle

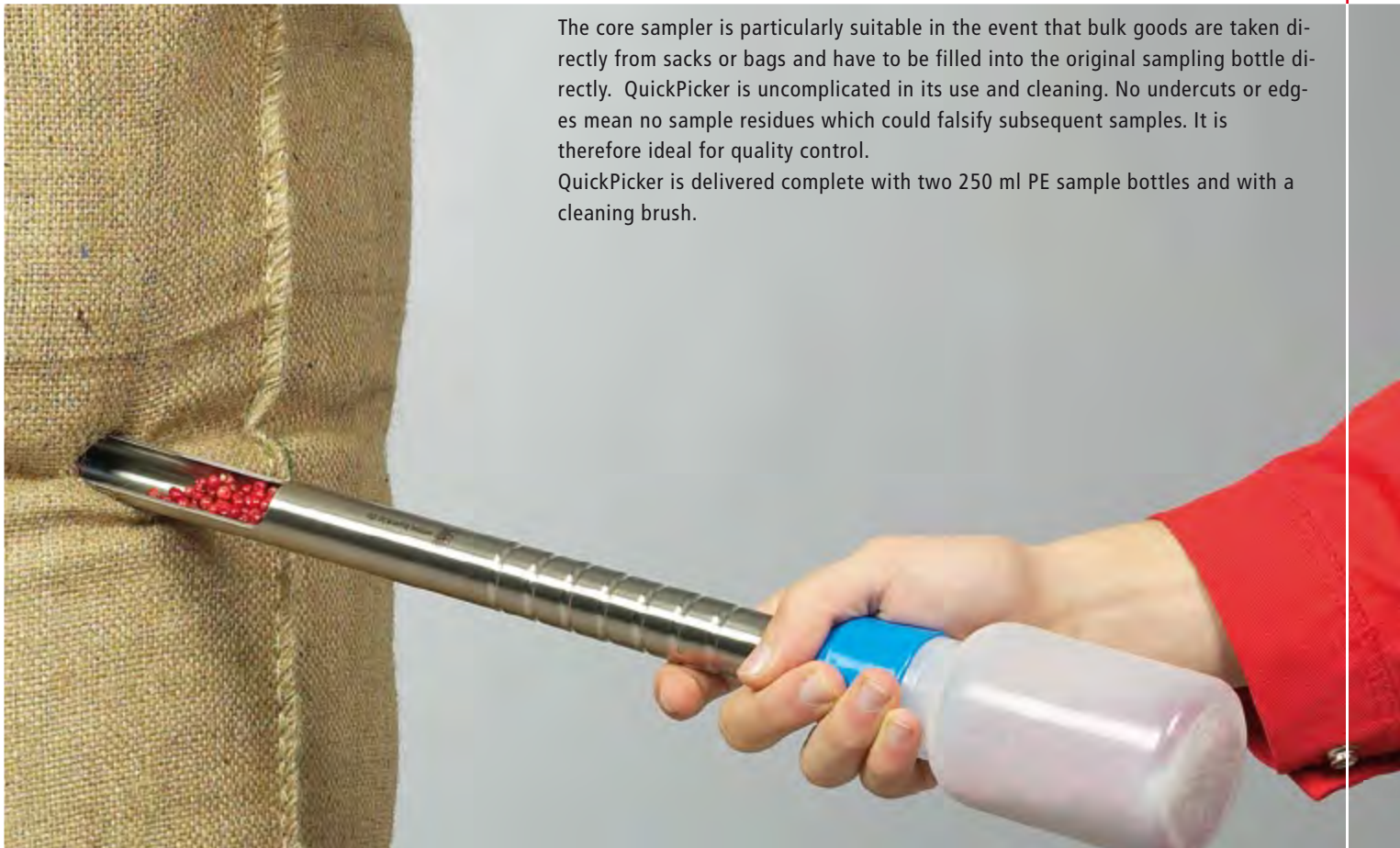

The core sampler is particularly suitable in the event that bulk goods are taken directly from sacks or bags and have to be filled into the original sampling bottle directly. QuickPicker is uncomplicated in its use and cleaning. No undercuts or edges mean no sample residues which could falsify subsequent samples. It is therefore ideal for quality control.

QuickPicker is delivered complete with two 250 ml PE sample bottles and with a cleaning brush.

QuickPicker | collects samples quickly and channels them directly into the sampling bottle

- ▶ Insertion depth 30 cm, diameter 25 mm
- ▶ Made of stainless steel (1.4404) approved for use with foodstuffs or PP transparent for aggressive chemicals, scratchproof, sterilisable, cost-effective
- ▶ Removable tip, easy cleaning
- ▶ Additional sampling bottles made of PE or PP available as accessories



QuickPicker V4A in use

i Corresponds to ISTA standard for seed lengths of 10 mm (wheat, oats) and 5 mm (flax, vetches).



close-it seals and close-it food lock seals ensure non-leak closure of the insertion point in bags and BigBags. See page 88.

QuickPicker

Material	Insertion depth mm	Total length cm	Volume ml	Price €	Quantity scale from	Scale price €	Item no.
V4A	300	50	75		5		5322-1300 N4
PP, transparent	300	50	75		5		5322-0300 N4

Accessories

Description	Price €	Item no.
Sampling bottle LDPE, 250 ml		0318-0250 N3
Sampling bottle PP, 250 ml		0319-0250 N3
Transport case for QuickPicker		3600-3000 N3
close-it control seals see page 88		